



Clinical Laboratories

**To:** *Clinical Laboratory staff*

**From:** *Laboratory Quality, Compliance and Safety*

**Date:** 9/27/2024

**Subject:** **Intelligent Locations User Interface Update Notice**

**Effective Date:** September 29, 2024


**Laboratory Section:** All Clinical laboratory section

**Summary:**

The Intelligent Locations Application will be permanently moving to a new user interface and the old will no longer be available starting on September 29, 2024 at 2:00am. Although most users have been utilizing the new interface already for several months now we wanted to provide a quick visual of where key functions are in the upgrade user interface. Although this update changes the view slightly all functions previously available and used remain as part of the application but are accessed in a slightly different location. Please refer to your applicable SOP UCMC 5956 which gives step by step instructions on how to use this interface.

**Old User Interface:**


- By selecting the “Get Details” icon circled in red below you can access an assets Details, Sensor History Data, Alarms, and Map View. This allowed you to review temperature reading, comment on alarms, disable alarms, and place a device in maintenance mode for extend service repairs.

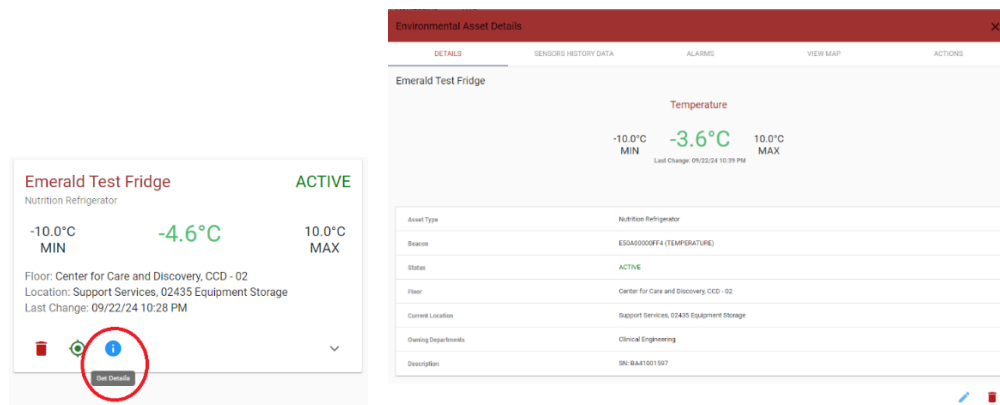



New User Interface:

- By selecting the “View Details” symbol circled below in orange you can access the details of an asset.
- By selecting the current location bubbles below a map of the location will popup.
- By selecting the three dots icon circled in red below you can access an assets, Action, Alarm, and Sensor History. Along with Configure Alarms and Turn Maintenance Mode On for extend service repairs.

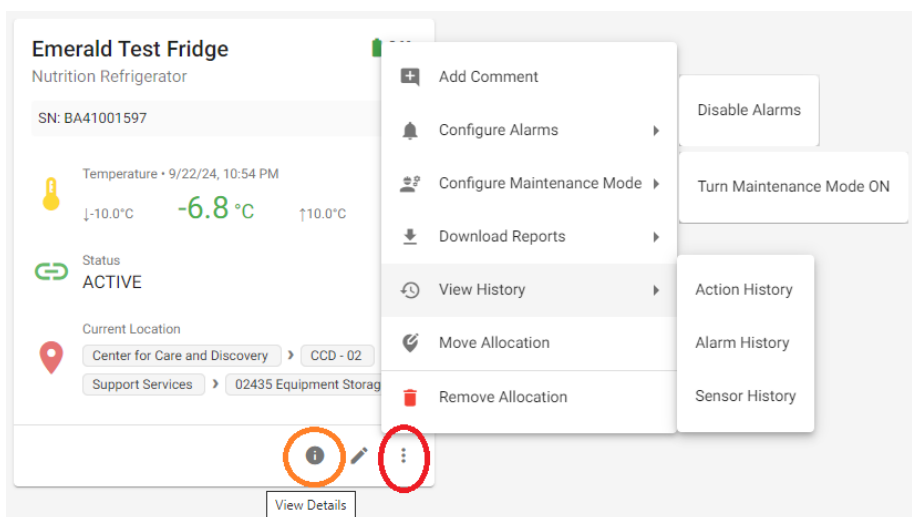

**Questions:** If there are any questions regarding the change, please contact: Lorena Chesebro at [Lorena.chesebro@uchicagomedicine.org](mailto:Lorena.chesebro@uchicagomedicine.org) or Marcus Holme at [marcus.holme.@uchicagomedicine.org](mailto:marcus.holme.@uchicagomedicine.org).